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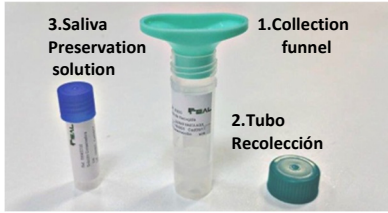
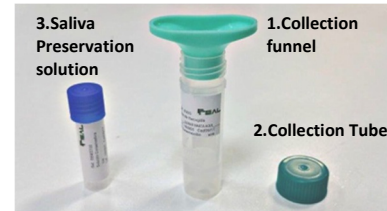
REAL SALIVA DNA Sample Collection Kit

Ref: RBMSAL01/100/500/1000

INTRODUCTION

This product is designed for simple and non-invasive collection and preservation saliva samples at room temperature. The saliva collection tube (2) is sent to the laboratory for DNA Isolation and analysis using the **REALPURE SALIVA KIT** or other purification kits.

The saliva preserved samples is stable for more than 2 years at room temperature.

PRESERVATION: Preservation between 15-25°C up to 3 years is recommended.

WARNINGS: Do not ingest the saliva preservation solution.

Keep away from children.

SAMPLE COLLECTION AND PRESERVATION PROCEDURES

1. Do not eat, chew gum, drink or smoke for 30 minutes prior to saliva collection.

2. Rinse your mouth with water and wait for 10 minutes.

3. Rub your cheeks and wiggle your tongue to stimulate the saliva secretion.

4. Place the collection funnel (1) in the saliva collection tube (2). Spit into the collection funnel until the liquid saliva (not including bubbles) in the collection tube reaches 2ml line.
If you notice excess bubbles eliminate them by gently tapping the tube on a hard surface.

5. Remove the tip of the saliva preservation solution (3).

6. Add the saliva preservation solution (3) to the collection tube.

7. Screw on the cap of the collection tube and close it tightly.

8. Shake well for 15 seconds to mix.

9. The saliva sample is now preserved and ready for storage, shipping or processing.