REAL DNA REMOVE PAQ 2X100ML

Ref.:RD055 (2 X 100 ML)

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1.INTRODUCTION

Nucleic acid amplification techniques have been a major breakthrough in the field of Molecular Biology: using methods as simple as the Polymerase Chain Reaction (PCR). A single DNA or RNA molecule can be rapidly detected from a wide variety of samples. However, on certain occasions and due to the extreme sensitivity of this technique, exogenous DNAs, plasmids or, above all, products of previous amplifications, can become the target DNA of subsequent amplification reactions, giving rise to false positives. To avoid contamination, there are control measures based on the use of enzymes such as Uracil DNA glycosylase, which involves modifications to the standard PCR method and an increase in the cost of the technique. Many other strategies have been suggested to minimise contamination, such as using filtered pipette tips, working in laminar flow cabinets, dividing the laboratory into pre-amplification and post-amplification areas, autoclaving the work material, etc. Some of these measures are not effective and others are unfeasible for many laboratories. DNA Remove is a detergent mixture specially formulated to remove DNA or RNA from work surfaces and laboratory materials in a very simple and effective way.

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3. GENERAL PROCEDURE

1. Spray the work surfaces and lab material with DNA Remove (you can also use a cloth wet in DNA Remove).
2. Wait for at least 3 minutes and wash with distilled water or ethanol.
3. To decontaminate the pipettes, submerge the autoclavable part in DNA Remove and wash with distilled water

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- DNA Remove is compatible with materials such as glass, PVC, PE and INOX. Avoid long time contact with surfaces like iron, nickel or copper.
- Use glasses and gloves to prevent splashes.
- In case of contact in eyes or skin wash with water.
- Store at room temperature. Avoid exposure to high temperature and intense light